

Water Sample Collection System

2000 mL water sampler for discreet and composite collections.

Construction

Ultra-high-molecular-weight polyethylene (UHMW), PVC, & stainless steel hardware.

1. Rope attachment point
2. Spring action opener
3. Handle
4. Seated plugs
5. Main body & lid (UHMW)
6. Isolated stainless steel weighted base

Sampler Dimensions

Main Body Including Lid
17.0" H, 4.5" Dia. (43.2 cm H, 11.5 cm Dia.)

Internal Sample Containment
10.75" H, 4.0" Dia. (27.3 cm H, 10.2 cm Dia.)

Shipping Dimensions & Weight

6" X 6" X 32" (15.2 cm X 15.2 cm X 81.3 cm)

13 pounds (5.9 kg)

Discreet Sample Deployment

Insert seated plugs and lower to desired depth. To open the sampler, give the attachment rope a good tug to engage the spring action. The seated plugs will pop out and sampling will commence.

Composite Sample Deployment

From surface: Remove seated plugs and lower through water column.

From given point in the water column: Deploy as if a discreet sample. Once at the desired starting depth, open sampler and then begin to lower through water column.

Pricing (USD, shipping included): \$2,160

Pricing is for U.S. domestic only (lower 48 States) and is subject to change without notice. For international inquiries, please contact us to discuss further.

To order, please enter in information below, sign, and email to watercare@ixom.com or contact us at **+1 866-437-8076**.

Company: _____

Name: _____

Phone: _____

Quantity: _____

Signature & Date: _____

